

## **CALIBREX DISPENSER**

### **520 Series (2,5,10ml)**



### **“HELPFUL CLEANING & DOSING HINTS”**

- Dispensers should be cleaned daily by rinsing in hot water.
- A more thorough clean should be undertaken weekly.  
(See cleaning protocol for detailed instructions)
- If disassembling the unit for cleaning, take care not to lose the small glass aspiration ball contained within the ceramic white disc.
- Should you lose the glass ball, replacements can be ordered from Interpath Services
- Components should be rinsed and soaked in hot water with a gentle detergent.
- When priming the unit for use after cleaning, dispense into a small clean container to ensure there are no air bubbles and recycle the solution back into the bottle.
- Should you need to send your unit back to Interpath Services for calibration or service, please ensure the Decontamination Certificate is included and correctly completed.
- Accurate dosing relies on consistent procedures particularly in speed and technique.
- A slow and steady dispensing technique ensures a complete uptake and delivery of liquid.
- Calibration checks should be conducted gravimetrically by weight using a digital balance according to ISO/AS 2162.2 protocol.
- Using glass measuring cylinders to calibrate dispensers should be avoided as their accuracy can vary.